

---

**User Guide**  
Elite-Cam  
Connection & Setup



## Elite-Cam Integrated USB Digital Camera

The Elite-Cam is a permanently-engaged USB digital camera solution for integration with Vision Engineering products.

### Health & Safety



Vision Engineering and its products conform to the requirements of the EC Directives on Waste Electrical and Electronic Equipment (WEEE) and Restriction of Hazardous Substances (RoHS).



---

## INTRODUCTION

---

Elite-Cam Integrated USB Digital Camera	i
Health & Safety	i

---

## CONTENTS

---

Pack contents	1
---------------	---

---

## GETTING STARTED

---

Assembly	2
Operation	2
General care and maintenance	2


---

## WARRANTY

---


### Pack contents

- 1 Elite-Cam viewing head
- 2 Software CD
- 3 Interconnect cable
- 4 Elite-Cam AMCap user guide (not shown)
- 5 Mantis Elite user guide (not shown)
- 6 Mantis Elite-Cam user guide (not shown)

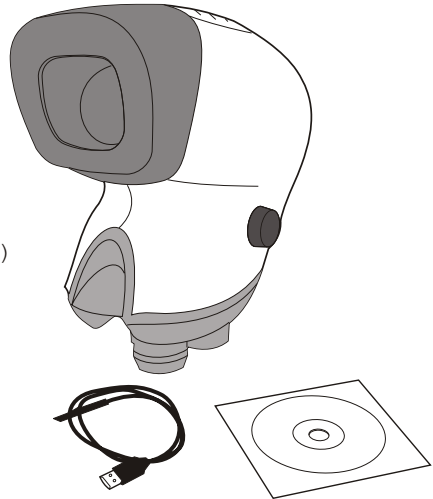

## Assembly

- ▶ Follow the Mantis Elite user guide to mount the viewing head **1** to the stand.
- ▶ Install software contained on the software CD **2**, as per Elite-Cam AMCap software user guide.
- ▶ Connect cable **3** to the viewing head **1** and then the chosen computer as per Elite-Cam AMCap software user guide.

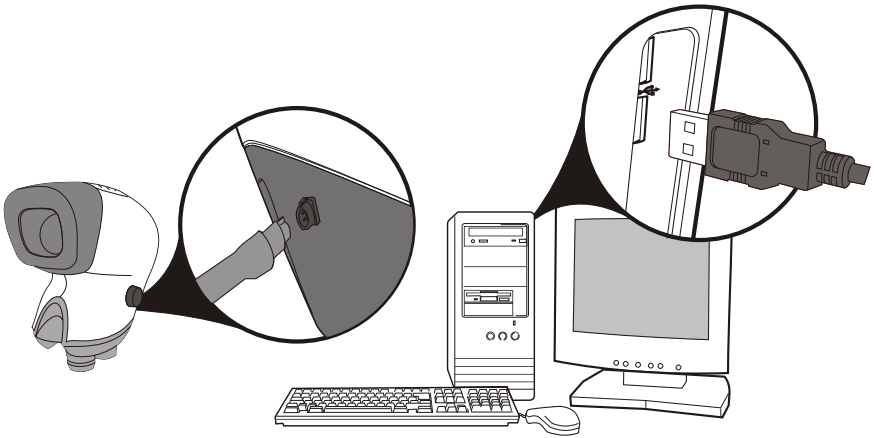

## Operation

- ▶ Refer to Mantis Elite user guide provided for general operation.
- ▶ Refer to Elite-Cam AMCap user guide to take an image or video capture of the subject.

## General care and maintenance

For all care and maintenance guidelines see the Mantis Elite user guide.



## **WARRANTY**

This product is warranted to be free from defects in material and workmanship for a period of one year from the date of invoice to the original purchaser.

If during the warranty period the product is found to be defective, it will be repaired or replaced at facilities of Vision Engineering or elsewhere, all at the option of Vision Engineering. However, Vision Engineering reserves the right to refund the purchase price if it is unable to provide replacement, and repair is not commercially practicable or cannot be timely made. Parts not of Vision Engineering manufacture carry only the warranty of their manufacturer. Expendable components such as fuses carry no warranty.

This warranty does not cover damage in transit, damage caused by misuse, neglect, or carelessness, or damage resulting from either improper servicing or modification by other than Vision Engineering approved service personnel. Further, this warranty does not cover any routine maintenance work on the product described in the user guide or any minor maintenance work which is reasonably expected to be performed by the purchaser.

No responsibility is assumed for unsatisfactory operating performance due to environmental conditions such as humidity, dust, corrosive chemicals, deposition of oil or other foreign matter, spillage, or other conditions beyond the control of Vision Engineering.

Except as stated herein, Vision Engineering makes no other warranties, express or implied by law, whether for resale, fitness for a particular purpose or otherwise. Further, Vision Engineering shall not under any circumstances be liable for incidental, consequential or other damages.

**For more information...**

Vision Engineering has a network of offices and technical distributors around the world. For more information, please contact your Vision Engineering branch, local authorised distributor, or visit our website.

Distributor

©

Vision Engineering Ltd.  
**(Manufacturing)**  
 Sand Road, Sand,  
 Woking, Surrey GU24 7ER, England  
 tel: +44 (0) 1483 248300  
 Email: generalinfo@visioneng.com

Vision Engineering Ltd.  
**(Commercial)**  
 Monument House, Monument Way West,  
 Woking, Surrey, GU21 5EN, England  
 tel: +44 (0) 1483 248300  
 Email: generalinfo@visioneng.com

Vision Engineering Inc.  
**(Manufacturing & Commercial)**  
 570 Danbury Road,  
 New Milford, CT 06776, USA  
 tel: +1 860 353 9776  
 Email: info@visioneng.com

Vision Engineering Inc.  
**(West Coast Commercial)**  
 745 West 14th Avenue,  
 Orange, CA 92665, USA  
 tel: +1 (714) 974 6966  
 Email: info@visioneng.com

Vision Engineering Ltd.  
**(Central Europe)**  
 Anton-Pendele-Str. 3,  
 82275 Emming, Deutschland  
 tel: +49 (0) 8141 40167-0  
 Email: info@visioneng.de

Vision Engineering Ltd.  
**(France)**  
 ZAC de la Nambiale, Av. de la Nambiale  
 91220 Le Plessis Palé, France  
 tel: +33 (0) 160 76 40 00  
 Email: info@visioneng.fr

Vision Engineering Ltd.  
**(Italy)**  
 Via Cesare Cantù, 9  
 20092 Cinisello Balsamo MI, Italia  
 tel: +39 02 6129 3518  
 Email: info@visioneng.it

Vision Engineering  
**(Brazil)**  
 tel: +55 11 4063 2206  
 Email: info@visioneng.com.br

Nippon Vision Engineering  
**(Japan)**  
 272-2 Saeba-cho, Tsuda-ku,  
 Yokohama-shi, 224-0054, Japan  
 tel: +81 (0) 45 935 1117  
 Email: info@visioneng.jp

Vision Engineering  
**(China)**  
 111, International Ocean Building,  
 720 Pudong Avenue, Shanghai,  
 200120, P.R. China  
 tel: +86 (0) 21 5036 7556  
 Email: info@visioneng.com.cn

Vision Engineering  
**(S.E. Asia)**  
 tel: +603 80700908  
 Email: info@visioneng.asia

Vision Engineering  
**(India)**  
 Email: info@visioneng.co.in

Visit our website:

# www.visioneng.com